

Effect on Lung Permeation of Various Formulation Efforts Invested in the Inhaled Antibiotic Apramycin as Measured in the *Ex Vivo* Isolated, Perfused, and Ventilated Rat Lung

Per Gerde

2025-12-10

Background

- The Isolated, ventilated and Perfused Lung of the rat (IPL), represents one versatile *ex vivo* model for use in preclinical drug development
- Integrated into the PreciseInhale[®] exposure platform, different aerosol formulations can be evaluated for lung PK and ventilation mechanics with a minimum of translational errors
- Four different formulations of the antibiotic Apramycin (Apr) was evaluated and compared for lung PK- and mechanics in the IPL system

APRINHA - New Inhalation Formulations for Improved Delivery of Antimicrobials



Duration: 36 month (Jan 2023 – Dec 2025)

Budget: 1 542 323 €

Project Objectives

Identify, characterize, and optimize a nanoparticle inhalation technology best suited for respiratory delivery of apramycin and other antimicrobials.

<https://www.jpiamr.eu/projects/aprinha/>

APRINHA - a European multi-center study

The consortium




Apramycin

Aminoglycoside antibiotic

Repurposed veterinary drug substance

Apramycin unique structure make it less susceptible to common aminoglycoside resistance mechanisms

DOI: 10.1021/jacs.9b11601

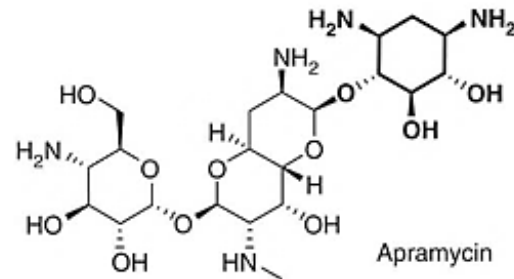


Physicochemical properties of Apramycin (free base)

Property	Value
MW	539.58 g/mole
Crystallinity	Crystalline
Melting point (Peak)	262 - 264 °C
pKa ¹	Basic: 7.7; 8.2; 8.4; 9.0; 9.6
Solubility in water ²	> 200 mg/mL below pH 8

¹Predicted values Marvin sketch 5.12.1

²A commercial veterinarian product at 200 mg/mL is available in some European countries. Experimental work performed at RISE indicates a solubility of at least 500 mg/mL.

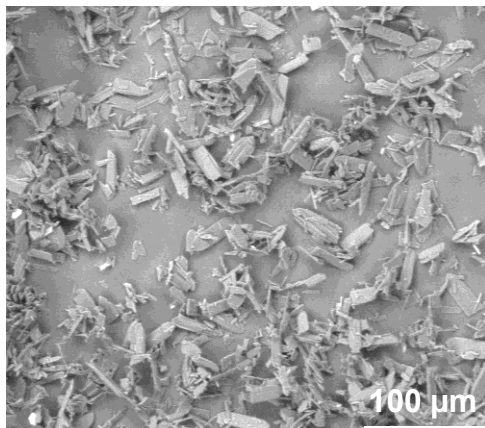

Inhalation formulation strategies for apramycin

Formulation 1 Base line reference Ref-Apr

- Apramycin solution (75 mg/mL, pH 5.5-6.0) for comparison of formulation performance

Solution

RI
SE

Formulation 2 Dextran-based NP KuDa-Apr

- Controlled release
- Biofilm penetration

Suspension

bio
cidetec>

Formulation 3 Micronized apramycin DSPC-Apr

- Decreased dissolution rate
- Low excipient to API

Powder


RI
SE

Formulation 4 Apramycin: Ga complex Ga-Apr



- Increased lung retention time
- Nanoparticle

Powder

Inserm

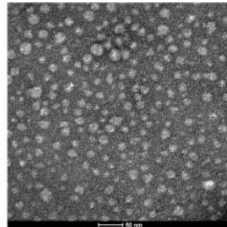

TEM. Dn = 13 ± 3 nm

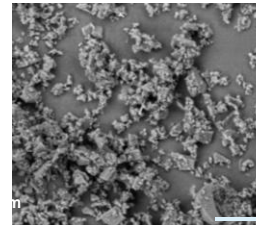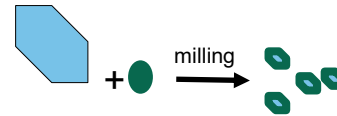

The PreciseInhale® Inhalation Switchboard

Aerosol sources

Dry powder

Clinical inhalers

Nebulized solutions

PreciseInhale®



Exposure modules

DissolvIt® – *in vitro*

- Non-biological dissolution and absorption testing
- Physiology-like conditions with single pass perfusion
- $C_{\max}$ and $T_{\max}$ PK profile curves

XposeALI® – *in vitro*

- Lung-like exposures of lung cells in culture
- Permeability assessment

Isolated Perfused Lung (IPL) – *ex vivo*

- Lung absorption rate and retention data
- <15% SD in variability during repetitive dosing
- High-resolution data on lung PK- and mechanics


Intratracheal inhalation in rodents – *in vivo*

- Direct exposure lungs avoiding GI exposure
- Spontaneous breathing animals – physiological data

Nose-Only inhalation in rodents – *in vivo*

- Precision by design – one animal at the time
- Individual control of exposure data

Isolated, ventilated and perfused lung of the rat (IPL)

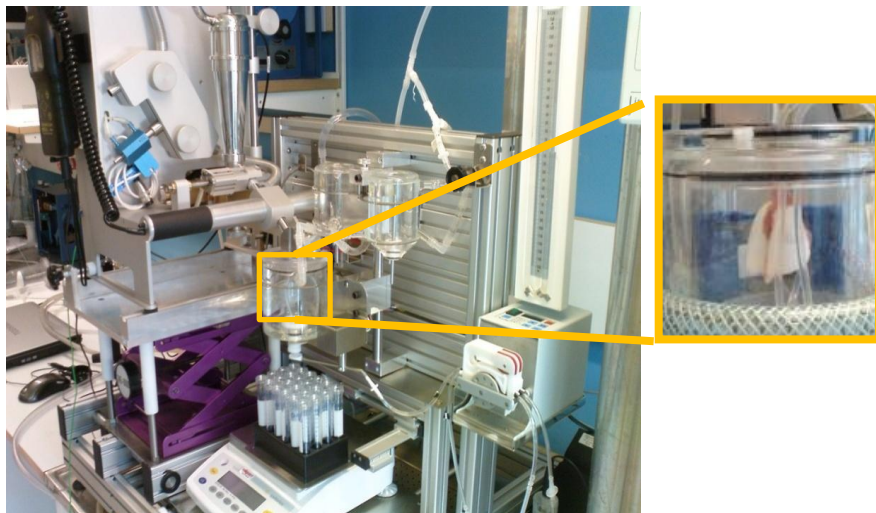


- Negative-pressure ventilated rat lung
- Aerosol exposure via a tracheal catheter
- Single-pass perfusion
- Two hour perfusate sampling time
- Monitoring of lung mechanics
- Over-all mass balance for dosing accuracy

Consistent Exposures with all Four Formulations of Apramycin


Formulation	MMAD (μm)	GSD	Deposition (μg , $\pm\text{SD}$)	Fraction of inhaled deposited (MPPD) Normalized regional deposition		
				Total of inhaled	Bronchial	Alveolar
1. Ref-Apr Solution	2.8	1.9	133 $\pm$ 9%	0.42	0.41	0.59
2. KuDa-Apr Suspension	3.6	2.2	151 $\pm$ 15%	0.49	0.51	0.49
3. DSPC-Apr Powder	3.0	2.2	117 $\pm$ 11%	0.43	0.46	0.54
4. Ga-Apr Powder	3.0	1.8	144 $\pm$ 12%	0.45	0.43	0.57

Distinctive Differences in Lung PK of Tested Formulations

A. Dose-normalized concentration in perfusate

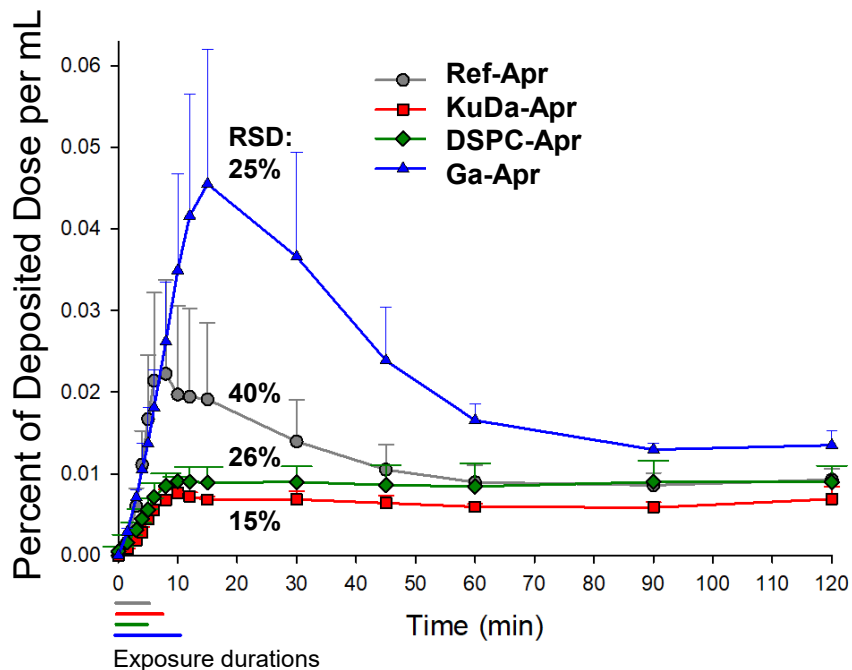

B. Fraction of deposited retained in lung

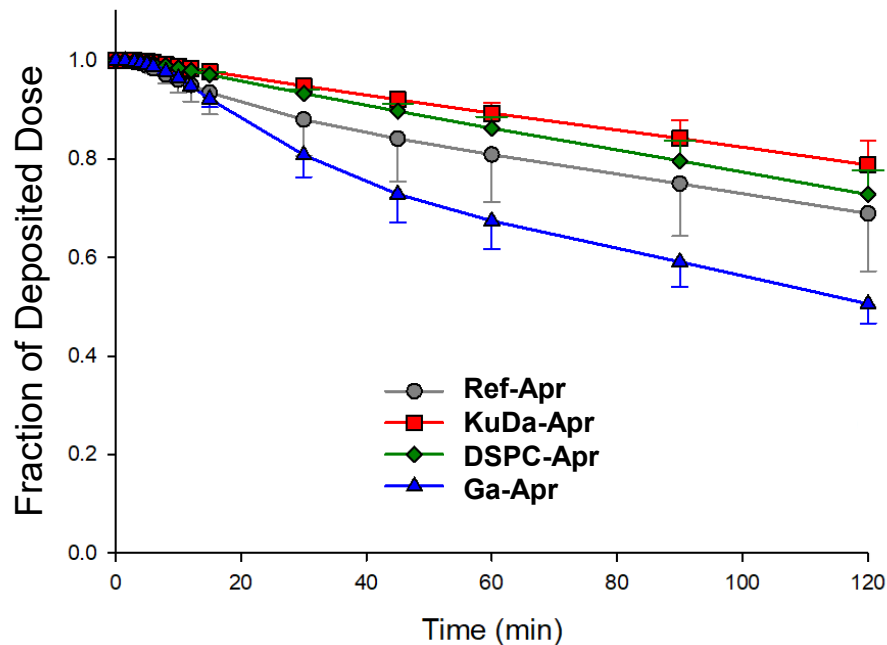

Key Results of Lung PK for Studied Formulations

Formulation	C _{max} (ng/mL)	T _{max} (min)	t _{1/2} (h)	Fraction retained at 2h	Fractional distribution at 2h post exposure			
					Lung	ELF	Perfusate	
1. Ref-Apr Solution	33 ± 50%	6.7 ± 17%	4.1 ± 47%	0.69 ± 17%	0.56 ± 19%	0.13 ± 20%	0.31 ± 38%	Ref
2. KuDa-Apr Suspension	13 ± 11%	9.3 ± 12%	6.0 ± 23%	0.79 ± 6%	0.55 ± 6%	0.24 ± 8%	0.21 ± 23%	Slower
3. DSPC-Apr Powder	21 ± 20%	7.0 ± 62%	4.5 ± 24%	0.73 ± 7%	0.55 ± 11%	0.14 ± 17%	0.31 ± 22%	Slower
4. Ga-Apr Powder	71 ± 47%	14 ± 12%	1.9 ± 14%	0.51 ± 8%	0.34 ± 15%	0.17 ± 35%	0.49 ± 8%	Faster

- Compared to Ref-Apr, KuDa-Apr and DSPC-Apr had an expected slower permeation in the lungs
- In contrast, Ga-Apr permeated the rat lung markedly faster than Ref-Apr


The Pulmonary Absorption Process of a iBCS Class III Drug

The Major Steps:

I Dissolution / Desorption



II Permeation

III Perfusion

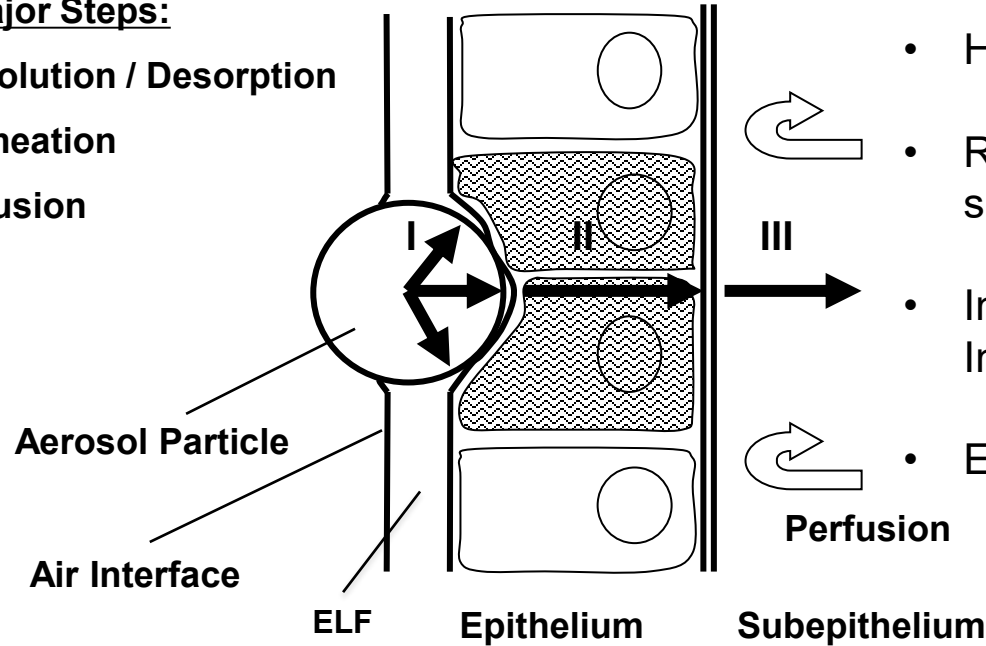

- High solubility, low permeability
- Rate control in step II; step I can only slow overall permeation
- Increase seen in Ga-Apr permeation! Increased barrier permeability?
- Evidence?

Effect of Apramycin Formulation Exposures on Lung Ventilation and Perfusion Rates

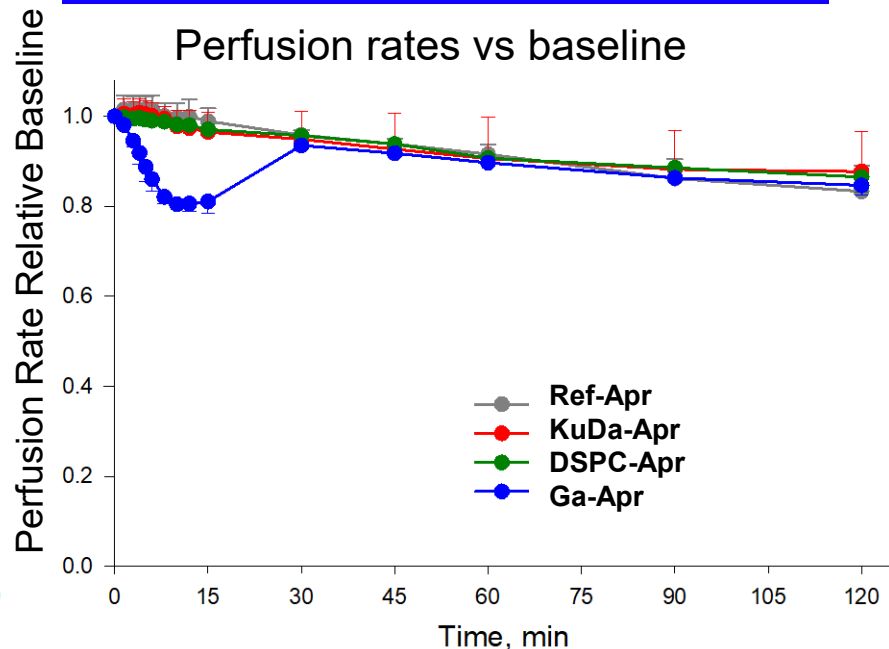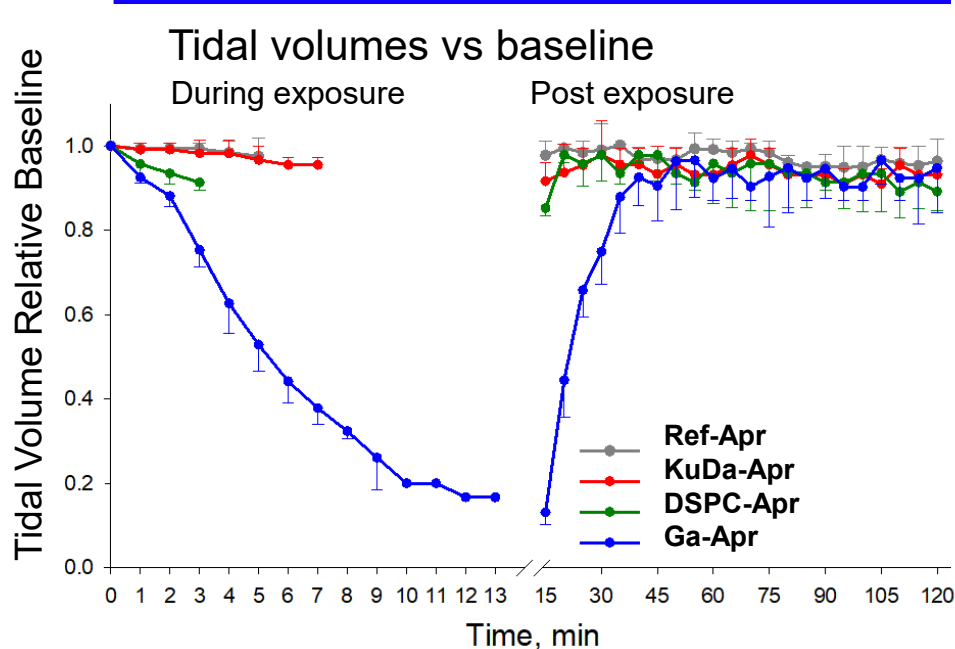

- Ref-Apr, KuDa-Apr, and DSPC-Apr exposures had no noticeable effect on lung physiology
- Ga-Apr caused a transient decrease in tidal volumes and perfusion rates for about 30-45 min
- No observable effect on lung physiology at the end of experiment

Conclusions

- For the iBCS Class III drug apramycin (high solubility and low permeability), it is possible to alter lung PK by using various engineered formulations
- Two formulations, KuDa-Apr and DPSC-Apr, slowed overall lung permeation compared to Ref-Apr
- One formulation, Ga-Apr, increased lung permeation compared to Ref-Apr, most likely by affecting the barrier function of the lung itself
- Low variability dosing with the PreciseInhale[®] system allow minimum number of animals to be used (n=3), still providing mechanistic understanding of lung PK (Reduce, Refine)

Thank you!

Inhalation Sciences

Ewa Selg

RISE

Sofie ALvér

Denny Mahlin

Bo Lassen

Martin Bolin

Mehrnaz Shaali

Marcus Sjödin

Puneet Rinwa

Poitiers, INSERM

Frédéric Tewes

Anissa Benkhaled

Vincent Aranzana-Climent

Sandrine Marchand

CIDETEC, BIO GIPUZKOA

Julie Movellan

José Carloz Ibañez

Damien Dupin

Iraida Loinaz

Fraunhofer, ITEM

Dorothee Winterberg

Vanessa Neuhaus

Katharina Schwarz

Karolin Sandra Webs

LIOS

Edgars Liepins